

WHITEPAPER

(READY-TO-PREP) RTP® Pathogen Kit for manual DNA/RNA extraction from pathogens

One of the greatest public health challenges of the 21st century are infectious diseases, due to complex and interrelated factors such as population growth, climate change, and geopolitics. Accurate diagnostic tools are needed for early detection of infectious diseases to prevent disease outbreaks. These diagnostic tools provide vital information for epidemiological models, disease tracking, treatment selection, and epidemic prevention. Diagnosis of infectious diseases at the genomic level using nucleic acid biomarkers has proven to be the most effective approach to date.

A nucleic acid extraction method specific to the sample material is of great importance for the success and informative value of sensitive downstream methods, such as amplification techniques. In human diagnostics, a wide variety of sample matrices are tested for pathogens and infectious diseases, e.g., swabs, tissue, stool samples or body fluids. Following sample extraction, various PCR techniques are used to detect a variety of pathogens of viral and bacterial origin. A particularly popular tool for straightforward sample analysis is multiplex PCR, which can detect viral DNA and RNA as well as bacterial DNA in a single assay. This method is frequently used for the diagnosis of respiratory infections. For the diverse sample matrices of human diagnostic and sophisticated downstream assays, the RTP® Pathogen Kit is the ideal application for sample extraction. Different nucleic acid types of viral and bacterial origin are extracted simultaneously. In addition, the RTP® Pathogen Kit offers a unique 1-step lysis, which increases reproducibility and helps to minimize pipetting steps in manual sample extraction.

RTP® - Ready-To-Prep

The RTP® Pathogen Kit greatly simplifies DNA/RNA purification from pathogen-containing samples. Everything needed for DNA/RNA sample preparation - lysis buffer, Proteinase K, lysozyme and carrier nucleic acids - is provided in a lyophilized format in an Extraction Tube (Fig. 1). The lyophilized buffer is ready-to-use and stable at room temperature. The kit is CE-IVD certified and complies with REGULATION (EU) 2017/746 on in vitro diagnostic medical devices.

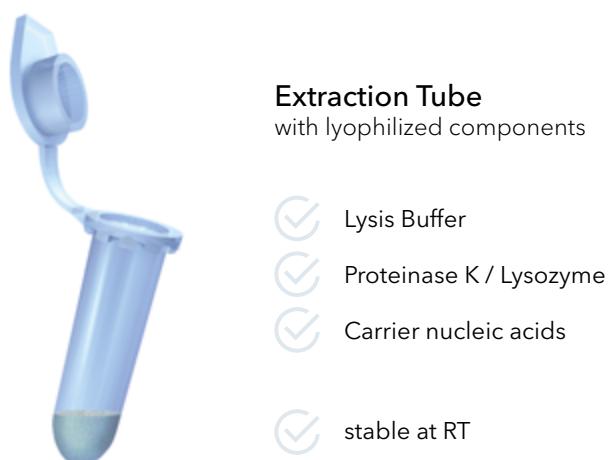


Figure 1: The Extraction Tube of the RTP® Pathogen Kit contains all components for sample lysis in a lyophilized buffer.

WHITEPAPER

EXTRACTION PROCEDURE

The one-step sample lysis takes place under non-chaotropic conditions, this gentle sample treatment allows potential inhibitors in the sample to be removed while nucleic acids remain intact. Due to the carrier nucleic acids in the extraction buffer, the recovery of small amounts of target sequences is enhanced.

After heat induced lysis, samples are transferred to a spin column-based procedure with a total preparation time of only 20 minutes. Impurities are removed by repeated washing steps, finally the purified nucleic acids can be eluted in a small volume of elution buffer.



Fig. 2: Invitek extraction kit

FAST AND CONVENIENT HANDLING

The RTP® technology reduces hands-on time and plastic waste while increasing reproducibility and safety. Up to 60% fewer pipetting steps and pipet tip consumption can be achieved with the one-step lysis compared to conventional kits (Fig. 3). Due to the broad-spectrum specification of the RTP® Pathogen Kit in terms of sample material and target nucleic acid, the kit can be used ubiquitously in the laboratory and enables a wide range of applications.

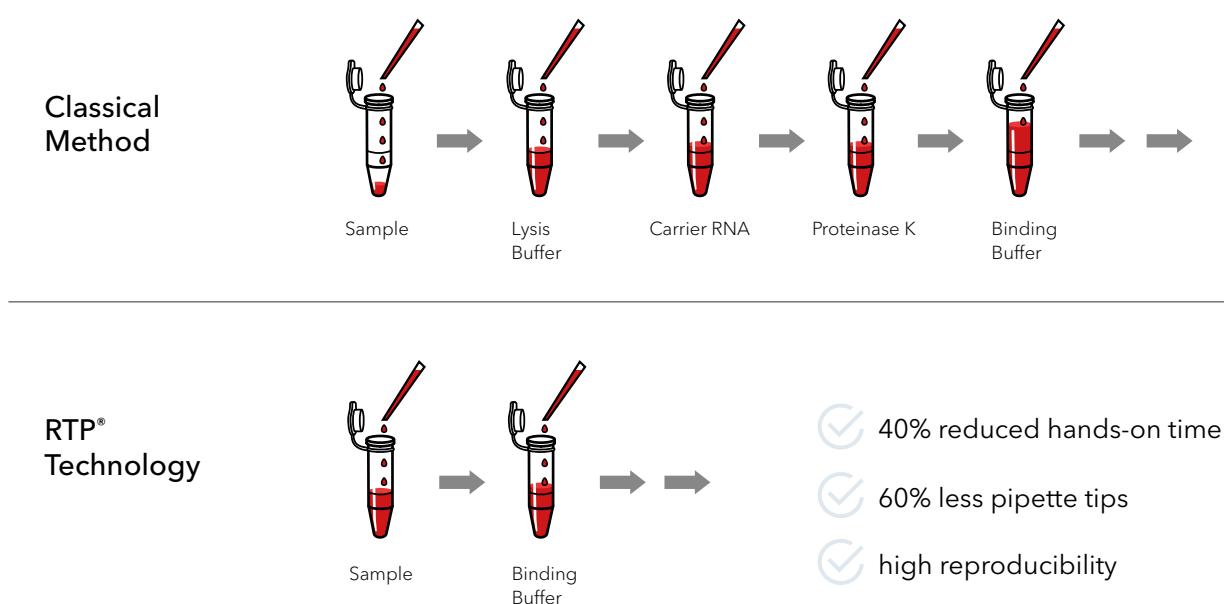


Figure 3: Comparison of a classical spin column extraction method with the RTP® technology. Due to a 1-step lysis, hands-on time and pipetting steps are reduced significantly during lysis while ensuring highest reproducibility.

Applications using the RTP® kit

For many years, Invitek's products have been used for the detection of infectious diseases, e.g., respiratory RNA viruses, DNA viruses such as adenovirus or herpes and various bacterial species. Due to the high-quality nucleic acids obtained with the extraction protocol classical and advanced molecular biological methods can be applied like multiplex qPCR and qRT-PCR and Next-Generation-Sequencing. Apart from diagnostic assays, the RTP® kit already has been used for diverse microbiome profiling studies.

WHITEPAPER

1. DNA extraction from bacteria

The literature encompasses diverse applications for isolating bacterial DNA, such as identifying human diseases, conducting genetic analyses of bacterial species, and detecting bacteria in food and environmental samples.

Table 1: Applications of the RTP® Pathogen Kit for bacterial detection in diagnostics and research

Pathogen	Starting material	Detection system
<i>E.coli</i> , <i>EHEC</i> , <i>EPEC</i>	Bacterial enrichment culture	Illumina whole-genome resequencing (GATC Biotech, Eurofins Genomics, Luxembourg) ¹ , GeneDisc® arrays (Pall, Port Washington, USA) ²
<i>Helicobacter pylori</i>	Bacterial enrichment culture	Pacific Biosciences RS sequencing technology (Pacific Biosciences, Menlo Park, USA) ³
<i>Legionella pneumophila</i>	Water, lung tissue or sputum	in-house PCR ^{4,5}
<i>Mycoplasma gallisepticum</i> , <i>Mycoplasma synoviae</i> , <i>Mycoplasma pneumoniae</i>	Bacterial culture, oropharyngeal swabs	in-house PCR ^{6,7,8}
<i>Yersinia enterocolitica</i>	Bacterial culture	in-house PCR ⁹
<i>Pseudomonas aeruginosa</i>	Sputum	in-house PCR ¹⁰
<i>Salmonella enterica</i>	Bacterial culture	in-house PCR, microarray, multilocus sequence typing ^{11,12}
<i>Vibrio cholerae</i> , <i>Vibrio vulnificus</i> , <i>Vibrio parahaemolyticus</i> , other <i>Vibriaceae</i>	Bacterial culture	in-house PCR ^{13,14,15,16,17}
<i>Treponema</i> spp.	Serum and CSF samples, brain tissue	in-house PCR ¹⁸
<i>MRSA</i>	Bacterial culture, bronchoalveolar lavage, throat swab/nasal swab	FTD-RP33 (Fast-Track Diagnostics, Luxembourg) ¹⁹ , EVIGENE MRSA detection kit (EVIGENE; Statens Serum Institut, Copenhagen, Denmark) ²⁰
<i>Bacillus cereus</i>	Bacterial culture	in-house PCR ²¹
<i>Mycobacterium tuberculosis</i>	Sputum	Genotype-MTBDR assay (Hain Life-science, Nehren, Germany) ²²
<i>Chlamydia trachomatis</i>	Urogenital swabs	in-house PCR ²³
<i>Periodontal bacteria</i> (<i>Porphyromonas gingivalis</i> , <i>Aggregatibacter actinomycetemcomitans</i> , <i>Tannerella forsythia</i> , <i>Treponema denticola</i> , <i>Prevotella intermedia</i>)	Paper points, saliva samples, plaque samples	in-house PCR ²⁴ , micro-IDent microorganism kit Hain Lifescience GmbH, Nehren, Germany ²⁵
<i>Environmental bacteria</i> (red sea bacterial strains)	Bacterial culture	in-house PCR ²⁶

Furthermore, the RTP® Pathogen Kit was used for microbiome analysis in human and plant samples:

Table 2: Applications of the RTP® Pathogen Kit for microbiome analysis

Target microbiome	Starting material	Downstream application
Gut	Bacterial culture	in-house PCR ^{27,28}
Skin	Skin swabs	in-house PCR ²⁹
Vaginal	Vaginal swabs	Next Generation Sequencing ³⁰
Oral	Paper points, saliva, oral biofilms from membrane filters	Pyrosequencing ³¹ , sequencing ³²
Plant	Potato, carrot, beet, neep, topinambur	in-house PCR ³³

WHITEPAPER

2. DNA/RNA extraction from viruses

Here, we present a listing of literature wherein viral RNA and DNA isolation is documented for the purpose of detecting viral diseases, including respiratory or tropical diseases.

Table 3: Applications of the RTP® Pathogen Kit for viral detection in diagnostics and research

Pathogen	Starting material	Detection system
Viral RNA		
SARS-CoV-2 and other Corona viruses	Nasopharyngeal swab, bronchoalveolar lavage	in-house PCR ^{34,35} , FTD-RP33, FTD-RP21 plus (Fast-Track Diagnostics, Luxembourg) ^{19,36,37,38}
Influenza A/B/C, H1N1, Parainfluenza virus	Nasopharyngeal swab, bronchoalveolar lavage	FTD-RP33, FTD-RP21 plus (Fast-Track Diagnostics, Luxembourg) ^{19,36,37,38} , RealStar® Influenza RT PCR Kit 1.0 (Altona Diagnostics, Hamburg, Germany) ³⁹
Respiratory syncytial virus (RSV)	Sputum, serum, nasal swab, throat swab	FTD-RP33, FTD-RP21 plus (Fast-Track Diagnostics, Luxembourg) ^{19,36,37,38} , RSV Serotype A and B kits (Roboscreen, Leipzig, Germany) ⁴⁰
Zika virus	Serum, urine samples	in-house PCR ⁴¹
Crimean-Congo hemorrhagic fever virus	Ticks	in-house PCR, RealStar CCHFV RT-PCR Kit 1.0 (Altona Diagnostics, Hamburg, Germany) ⁴²
Rotavirus	Virus culture	in-house PCR ⁴³
Usutu virus	Mosquitoes, tissue samples	in-house PCR ⁴⁴
Phlebovirus and Massilia virus	Sandflies	in-house PCR ⁴⁵
Tick-borne encephalitis virus (TBEV)	Tissue sample (spleen)	in-house PCR ⁴⁶
Hepatitis C, E	Liver tissue	INNO-LiPA™ HCV II kit (Innogenetics, Ghent, Belgium) ⁴⁷ , in-house PCR ⁴⁸
Viral DNA		
Adenovirus	Water, sediments, stool suspension, sputum, serum, urine, bronchoalveolar lavage, throat swab/nasal swab	FTD-RP33, FTD-RP21 plus (Fast-Track Diagnostics, Luxembourg) ^{19,36,37,38} , Adenoplex® multiplex PCR kit (Gen-Probe Prodesse Inc., Waukesha, USA) ⁴⁰
Human Cytomegalovirus (CMV), Epstein Barr Virus (EBV), Human Papilloma Virus (HPV)	Tissue	in-house PCR ^{49,50,51}
Herpes Simplex Virus 1 & 2, Varicella-zoster virus (VZV)	Cell culture supernatant, CSF, vesicle fluid samples	in-house PCR ^{52,53,54,55}
Hepatitis B	Oral fluid samples	in-house PCR, Abbott Real Time HBV (Abbott, Abbott Park, Illinois, USA) ⁵⁶
Equid herpesviruses	Cell culture supernatant, swabs	in-house PCR ⁵⁷
African swine fever virus	Insect tissue	in-house PCR ⁵⁸
BK and JC polyomaviruses	Tissue	in-house PCR ⁵⁹
Orthopoxvirus	Blotting paper	in-house PCR ⁶⁰

WHITEPAPER

RTP® Pathogen Kit in comparison to other target specific extraction kits

Previously, there were several RTP® kits specifically offered for bacteria or viruses (RTP® Bacteria DNA Mini Kit, RTP® DNA/RNA Virus Mini Kit, RTP® Mycobacteria Kit). For easier and more flexible handling, Invitek offers the RTP® Pathogen Kit with different protocols for bacteria and viruses combined in one kit.

To compare the performance of the RTP® Pathogen Kit and RTP® DNA/RNA Virus Mini Kit, RNA was prepared from 0.1 µl Influenza A virus stock spiked in 200 µl serum. Three replicates each were purified with both extraction kits. PCR was performed with SureFast® Influenza A Virus Kit (Congen, Berlin, Germany) according to the manufacturer's instructions. Results show (Fig.4) equivalent results for both extraction kits. Additionally, to compare the performance of the RTP® Pathogen Kit and RTP® Bacteria DNA Mini Kit, three replicates each were extracted from a *Bacillus subtilis* overnight culture. In the subsequent kit performance assay, PCR was done specifically for the internal control: ECD = Extraction Control DNA. Results show (Fig.5) equivalent results for both extraction kits.

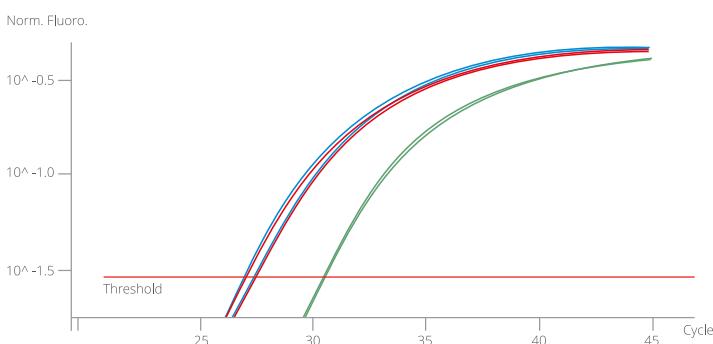


Figure 4: PCR assay SureFast® Influenza A (Congen, Berlin, Germany), FAM Channel. RNA was extracted from Influenza A spiked serum with RTP® Pathogen Kit and RTP® DNA/RNA Virus Mini Kit. Blue: extraction with RTP® Pathogen Kit. Red: extraction with RTP® DNA/RNA Virus Mini Kit. Green: PTC.

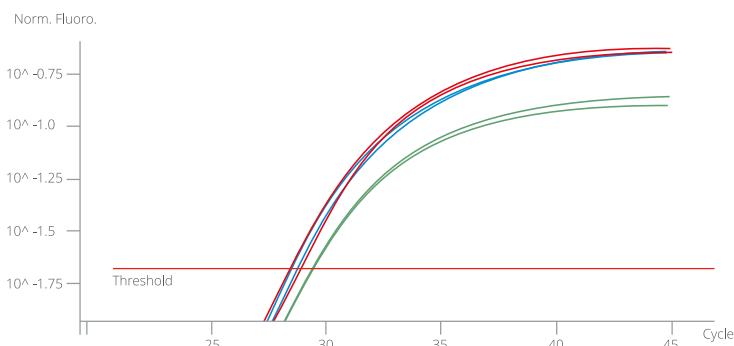


Figure 5: PCR assay specific for ECD, FAM Channel. Blue: extraction with RTP® Pathogen Kit. Red: extraction with RTP® Bacteria DNA Mini Kit. Green: PTC.

Conclusion

Invitek Diagnostics offers a reliable and convenient solution for nucleic acid extraction from bacteria, and viruses, which has been shown in many peer-reviewed publications.

- Isolation of viral DNA/RNA and bacterial DNA
- Suitable for all clinically relevant starting materials (e.g. serum, plasma, stool suspension, bacterial enrichment culture, tissue)
- Ready-to-Prep (RTP®) Technology: Pre-filled Extraction Tubes for one-step sample lysis, kits can be stored at room temperature
- Perfect for manual use: convenient handling and maximum reproducibility
- CE-IVD certified, recommended for in vitro diagnostic use

For automated use, we recommend the Invitek Universal Kit series which can be easily automated on several automated platforms, such as the KingFisher™ Flex.

WHITEPAPER

- 1 Stock culture heterogeneity rather than new mutational variation complicates short-term cell physiology studies of *Escherichia coli* K-12 MG1655 in continuous culture. Nahku R, Peebo K, Valgepea K, Barrick JE, Adamberg K, Vilu R. *Microbiology (Reading)*. 2011 Sep;157(Pt 9):2604-2610.
- 2 Virulence gene profiling of enterohemorrhagic (EHEC) and enteropathogenic (EPEC) *Escherichia coli* strains: a basis for molecular risk assessment of typical and atypical EPEC strains. Bugarel M, Martin A, Fach P, Beutin L. *BMC Microbiol*. 2011 Jun 21;11:142.
- 3 Comparing the genomes of *Helicobacter pylori* clinical strain UM032 and Mice-adapted derivatives. Khosravi Y, Rehvathy V, Wee WY, Wang S, Baybayan P, Singh S, Ashby M, Ong J, Amoyo AA, Wee SS, Choo SW, Perkins T, Chua EG, Tay A, Marshall BJ, Loke MF, Goh KL, Pettersson S, Vadivelu J. *Gut Pathog*. 2013 Aug 19;5(1):25.
- 4 Prevalence of sequence types among clinical and environmental isolates of *Legionella pneumophila* serogroup 1 in the United States from 1982 to 2012. Kozak-Muiznieks NA, Lucas CE, Brown E, Pondo T, Taylor TH Jr, Frace M, Miskowski D, Winchell JM. *J Clin Microbiol*. 2014 Jan;52(1):201-11.
- 5 Application of TaqMan low-density arrays for simultaneous detection of multiple respiratory pathogens. Kodani M, Yang G, Conklin LM, Travis TC, Whitney CG, Anderson LJ, Schrag SJ, Taylor TH Jr, Beall BW, Breiman RF, Feikin DR, Njenga MK, Mayer LW, Oberste MS, Tondella ML, Winchell JM, Lindstrom SL, Erdman DD, Fields BS. *J Clin Microbiol*. 2011 Jun;49(6):2175-82.
- 6 Mycoplasma gallisepticum and Mycoplasma synoviae express a cysteine protease CysP, which can cleave chicken IgG into Fab and Fc. Cizelj I, Bercic RL, Dusanic D, Narat M, Kos J, Dovc P, Bencina D. *Microbiology*. 2011 Feb;157(Pt 2):362-72.
- 7 Comparison of nucleic acid extraction methods for the detection of *Mycoplasma pneumoniae*. Thurman KA, Cowart KC, Winchell JM. 2009. *Diagn Microbiol Infect Dis* 65:435-438.
- 8 *Mycoplasma pneumoniae* from the Respiratory Tract and Beyond. Waites KB, Xiao L, Liu Y, Balish MF, Atkinson TP. *Clin Microbiol Rev*. 2017 Jul;30(3):747-809.
- 9 Characterization of a *Yersinia enterocolitica* biotype 1A strain harbouring an ail gene. Kraushaar B, Dieckmann R, Wittwer M, Knabner D, Konietzny A, Mäde D, Strauch E. *J Appl Microbiol*. 2011 Oct;111(4):997-1005.
- 10 Application of quantitative PCR to the diagnosis and monitoring of *Pseudomonas aeruginosa* colonization in 5-18-year-old cystic fibrosis patients. Typhaine Billard-Pomares, Stéphanie Herwegh, Nathalie Wizla-Derambure, Dominique Turck, René Courcol, and Marie-Odile Husson. *J. Med. Microbiol.*, Feb 2011; 60: 157 - 161.
- 11 Population structure of *Salmonella enterica* serovar 4,[5],12:b:- strains and likely sources of human infection. Toboldt A, Tietze E, Helmuth R, Junker E, Fruth A, Malorny B. *Appl Environ Microbiol*. 2013 Sep;79(17):5121-9.
- 12 Human infections attributable to the D-tartrate-fermenting variant of *Salmonella enterica* serovar Paratyphi B in Germany originate in reptiles and, on rare occasions, poultry. Toboldt A, Tietze E, Helmuth R, Fruth A, Junker E, Malorny B. *Appl Environ Microbiol*. 2012 Oct;78(20):7347-57.
- 13 Multiplex PCR for detection of virulence markers of *Vibrio vulnificus*. Bier N, Diescher S, Strauch E. *Lett Appl Microbiol*. 2015 May;60(5):414-20.
- 14 Virulence Profiles of *Vibrio vulnificus* in German Coastal Waters, a Comparison of North Sea and Baltic Sea Isolates. Bier N, Jäckel C, Dieckmann R, Brennholt N, Böer SI, Strauch E. *Int J Environ Res Public Health*. 2015 Dec 15;12(12):15943-59.
- 15 Genetic and phenotypic analysis of *Vibrio cholerae* non-O1, non-O139 isolated from German and Austrian patients. Schirm eister F, Dieckmann R, Bechlars S, Bier N, Faruque SM, Strauch E. *Eur J Clin Microbiol Infect Dis*. 2014 May;33(5):767-78.
- 16 Characterization of trh2 Harbouring *Vibrio parahaemolyticus* Strains Isolated in Germany. Bechlars S, Jäckel C, Diescher S, Wüstenhagen DA, Kubick S, Dieckmann R, Strauch E. *PLoS One*. 2015 Mar 23;10(3):e0118559.
- 17 Rapid identification and characterization of *Vibrio* species using whole-cell MALDI-TOF mass spectrometry. Dieckmann R, Strauch E, Alter T. *J Appl Microbiol*. 2010 Jul;109(1):199-211.
- 18 Increased occurrence of *Treponema* spp. and double-species infections in patients with Alzheimer's disease. Nemergut M, Bat kova T, Vigasova D, Bartos M, Hlozankova M, Schenk mayerova A, Liskova B, Sheardova K, Vyhalek M, Hort J, Laczo J, Kovacova I, Sitina M, Matej R, Jancalek R, Marek M, Damborsky J. *Sci Total Environ*. 2022 Oct 20;844:157114.
- 19 Spectrum of respiratory viral infections in liver disease patients with cirrhosis admitted in critical care unit. Bajpai V, Gupta E, Mitra LG, Kumar H, Maiwall R, Soni KD, Gupta A. *J Lab Physicians*. 2019 Oct-Dec;11(4):356-360.
- 20 Methicillin-resistant *Staphylococcus aureus* in cattle food chains - prevalence, diversity, and antimicrobial resistance in Germany. Tenhagen BA, Vossenkuhl B, Käsbohrer A, Alt K, Kraushaar B, Guerra B, Schroeter A, Fetsch A. *J Anim Sci*. 2014 Jun;92(6):2741-51.
- 21 Multiple-Locus Sequence Typing and Analysis of Toxin Genes in *Bacillus cereus* Food-Borne Isolates. Barbara Cardazzo, Enrico Negrisolo, Lisa Carraro, Leonardo Alberghini, Tomaso Patarnello, and Valerio Giaccone. *Appl. Envir. Microbiol.*, Feb 2008; 74: 850 - 860.
- 22 Performance of genotype-MTBDR test directly on clinical specimens. Gülden Yılmaz, Cafer Eroğlu, Halil Kurt, Melike Atase ver, Ahmet Saniç; *Journal of Microbiology and Infectious Diseases*; 2012; 2 (4): 135-141.

WHITEPAPER

- 23 Prevalence of Ureaplasma urealyticum, Mycoplasma hominis and Chlamydia trachomatis in symptomatic and asymptomatic patients. Cutoiu A, Boda D. *Biomed Rep.* 2023 Sep 1;19(4):74.
- 24 Oxidative Stress Parameters in Saliva and Its Association with Periodontal Disease and Types of Bacteria. Almerich-Silla JM, Montiel-Company JM, Pastor S, Serrano F, Puig-Silla M, Dasí F. *Dis Markers.* 2015;2015:653537.
- 25 Periodontopathogen profile of healthy and oral lichen planus patients with gingivitis or periodontitis. Ertugrul AS, Arslan U, Dursun R, Hakki SS. *Int J Oral Sci.* 2013 Jun;5(2):92-7.
- 26 Large scale patterns of antimicrofouling defenses in the hard coral *Pocillopora verrucosa* in an environmental gradient along the Saudi Arabian coast of the Red Sea. Wahl M, Al Sofyani A, Saha M, Kruse I, Lenz M, Sawall Y. *PLoS One.* 2014 Dec 8;9(12):e106573.
- 27 The Beta-Lactam Resistome Expressed by Aerobic and Anaerobic Bacteria Isolated from Human Feces of Healthy Donors. Vázquez-López R, Solano-Gálvez S, Álvarez-Hernández DA, Ascencio-Aragón JA, Gómez-Conde E, Piña-Leyva C, Lara-Lozano M, Guerrero-González T, González-Barrios JA. *Pharmaceuticals (Basel).* 2021 Jun 3;14(6):533.
- 28 Evaluation of inter-individual differences in gut bacterial isoflavone bioactivation in humans by PCR-based targeting of genes involved in equol formation. Braune A, Blaut M. *J Appl Microbiol.* 2018 Jan;124(1):220-231.
- 29 Degradation of benzo[a]pyrene by bacterial isolates from human skin. Sowada J, Schmalenberger A, Ebner I, Luch A, Tralau T. *FEMS Microbiol Ecol.* 2014 Apr;88(1):129-39.
- 30 Metagenomics Analysis Using Next Generation Sequencing of Vaginal Samples from Community Practices in the US. Budd WT, Harwich M, Bostwick GD, Meyers G, Dilts J, et al. (2015) *MOJ Cell Sci Rep* 2(2): 00020.
- 31 Stimulated and unstimulated saliva samples have significantly different bacterial profiles. Gomar-Vercher S, Simón-Soro A, Montiel-Company JM, Almerich-Silla JM, Mira A. *PLoS One.* 2018 Jun 1;13(6):e0198021.
- 32 Composition of microbial oral biofilms during maturation in young healthy adults. Langfeldt D, Neulinger SC, Heuer W, Staufenbiel I, Künzel S, Baines JF, Eberhard J, Schmitz RA. *PLoS One.* 2014 Feb 4;9(2):e87449.
- 33 Endophytic bacterial communities in peels and pulp of five root vegetables. Kõiv V, Arbo K, Maiväli Ü, Kisand V, Roosaare M, Remm M, Tenson T. *PLoS One.* 2019 Jan 11;14(1):e0210542.
- 34 SARS-CoV-2 and influenza virus coinfection among patients with severe acute respiratory infection during the first wave of COVID-19 pandemic in Bangladesh: a hospital-based descriptive study. Akhtar Z, Islam MA, Aleem MA, Mah-E-Muneer S, Ahmmmed MK, Ghosh PK, Rahman M, Rahman MZ, Sumiya MK, Rahman MM, Shirin T, Alamgir ASM, Banu S, Rahman M, Chowdhury F. *BMJ Open.* 2021 Nov 29;11(11):e053768.
- 35 Evaluation of recombinase-based isothermal amplification assays for point-of-need detection of SARS-CoV-2 in resource-limited settings. Ghosh P, Chowdhury R, Hossain ME, Hossain F, Miah M, Rashid MU, Baker J, Rahman MZ, Rahman M, Ma X, Duthie MS, Wahed AAE, Mondal D. *Int J Infect Dis.* 2021 Nov 7;114:105-111.
- 36 Respiratory Pathogens in Infants Diagnosed with Acute Lower Respiratory Tract Infection in a Tertiary Care Hospital of Western India Using Multiplex Real Time PCR. Sonawane AA, Shastri J, Bavdekar SB. *Indian J Pediatr.* 2019 May;86(5):433-438.
- 37 Acute Respiratory Illness in Rural Haiti. Kim YY, Lew JF, Keith B, Telisma T, Nelson EJ, Brantly AC, Chavannes S, Anilis G, Yang Y, Liu M, Alam MT, Rashid MH, Morris JG Jr, Madsen Beau De Rochars VE. *Int J Infect Dis.* 2019 Apr;81:176-183.
- 38 Acute respiratory viral infections in pediatric cancer patients undergoing chemotherapy. Benites EC, Cabrini DP, Silva AC, Silva JC, Catalan DT, Berezin EN, Cardoso MR, Passos SD. *J Pediatr (Rio J).* 2014 Jul-Aug;90(4):370-6.
- 39 Burden of influenza among hospitalized febrile children in Ghana. Hogan et al., *Influenza Other Respir Viruses.* 2017 Nov;11(6):497-501.
- 40 Detection of adenovirus and respiratory syncytial virus in patients with chronic obstructive pulmonary disease: Exacerbation versus stable condition. Kokturk N, Bozdayi G, Yilmaz S, Doğan B, Gulbahar O, Rota S, Tatlıcioğlu T. *Mol Med Rep.* 2015 Aug;12(2):3039-46.
- 41 Prolonged detection of Zika virus RNA in urine samples during the ongoing Zika virus epidemic in Brazil. de M Campos R, Cirne-Santos C, Meira GL, Santos LL, de Meneses MD, Friedrich J, Jansen S, Ribeiro MS, da Cruz IC, Schmidt-Chanasit J, Ferreira DF. *J Clin Virol.* 2016 Apr;77:69-70.
- 42 Crimean-Congo hemorrhagic fever virus clades V and VI (Europe 1 and 2) in ticks in Kosovo, 2012. Sherifi K, Cadar D, Muji S, Robaj A, Ahmeti S, Jakupi X, Emmerich P, Krüger A. *PLoS Negl Trop Dis.* 2014 Sep 25;8(9):e3168.
- 43 Identification of an avian group A rotavirus containing a novel VP4 gene with a close relationship to those of mammalian rotaviruses. Trojnar E, Sachsenröder J, Twardziok S, Reetz J, Otto PH, Johne R. *J Gen Virol.* 2013 Jan;94(Pt 1):136-142.
- 44 Putative Lineage of Novel African Usutu Virus, Central Europe. Cadar D, Bosch S, Jöst H, Börstler J, Garigliany MM, Becker N, Schmidt-Chanasit J. *Emerg Infect Dis.* 2015 Sep;21(9):1647-50.
- 45 Co-circulation of a novel phlebovirus and Massilia virus in sandflies, Portugal. Amaro F, Zé-Zé L, Alves MJ, Börstler J, Clos J, Lorenzen S, Becker SC, Schmidt-Chanasit J, Cadar D. *Virol J.* 2015 Oct 24;12:174.

WHITEPAPER

- 46 Flavivirus genome detected in Mus and Apodemus species sampled in illegal waste sites. Duh D, Varljen-Buzan E, Kunej U, Hasic S. Parasit Vectors. 2014 Apr 1;7(Suppl 1):P13.
- 47 Distribution of different hepatitis C virus genotypes in patients with hepatitis C virus infection. Bokharaei Salim F, Keyvani H, Amiri A, Jahanbakhsh Sefidi F, Shakeri R, Zamani F. World J Gastroenterol. 2010 Apr 28;16(16):2005-9.
- 48 Hepatitis E virus genotype 3 strains in domestic pigs. Cameroon. S de Paula V, Wiele M, Mbunkah AH, Daniel AM, Kingsley MT, Schmidt-Chanasit J. Emerg Infect Dis. 2013 Apr;19(4):666-8.
- 49 Cytomegalovirus and Langerhans Cell Histiocytosis: Is There a Link? Maliheh Khoddami, Seyed-Alireza Nadji, Paria Dehghanian, Mahsa Vahdatinia and Ahmad-Reza Shamshiri, Iran J Pediatr. 2016 April; 26(2):e673.
- 50 Detection of Epstein-Barr Virus DNA in Langerhans Cell Histiocytosis. Khoddami M, Nadji SA, Dehghanian P, Vahdatinia M, Shamshiri AR .Jundishapur J Microbiol. 2015 Dec 26;8(12):e2721.
- 51 Absence of Human Papillomavirus in Benign and Malignant Breast Tissue. Kazemi Aghdam M, Nadji SA, Alvandimanesh A, Maliheh M, Khademi Y. Iran J Pathol. 2019 Fall;14(4):279-283.
- 52 Comparison of different polymerase chain reaction methods for detection of herpes simplex virus types 1 and 2 encephalitis. Altuglu I, Zeytinoglu A, Sirin H, Yuceyar N, Erensoy S. Eur J Clin Microbiol Infect Dis. 2006 Oct;25(10):669-71.
- 53 A 12-year molecular survey of clinical herpes simplex virus type 2 isolates demonstrates the circulation of clade A and B strains in Germany. Schmidt-Chanasit J, Bialonski A, Heinemann P, Ulrich RG, Günther S, Rabenau HF, Doerr HW. J Clin Virol. 2010 Jul;48(3):208-11.
- 54 Novel varicella-zoster virus glycoprotein E gene mutations associated with genotypes A and D. Schmidt-Chanasit et al.; J. Clin. Microbiol., Jan. 2008, 46 (1); 325-327.
- 55 Molecular analysis of varicella-zoster virus strains circulating in Tanzania demonstrating the presence of genotype M1. Schmidt-Chanasit J, Olschläger S, Günther S, Jaeger G, Bleymehl K, Schäd SG, Heckel G, Ulrich RG, Doerr HW. J Clin Microbiol. 2008 Oct;46(10):3530-3.
- 56 Detection of occult hepatitis B in serum and oral fluid samples. Portilho MM, Nabuco LC, Villela-Nogueira CA, Brandão-Mello CE, Pilotto JH, Flores GL, Lewis-Ximenez LL, Lampe E, Villar LM. Mem Inst Oswaldo Cruz. 2018 Jan;113(1):62-65.
- 57 Detection of equid herpesviruses among different Arabian horse populations in Egypt. Azab W, Bedair S, Abdelgawad A, Eschke K, Farag GK, Abdel-Raheim A, Greenwood AD, Osterrieder N, Ali AAH. Vet Med Sci. 2019 Aug;5(3):361-371.
- 58 Trace amounts of African swine fever virus DNA detected in insects collected from an infected pig farm in Estonia. Herm R, Tummeleht L, Jürison M, Vilem A, Viltrop A. Vet Med Sci. 2020 Feb;6(1):100-104.
- 59 A globally applicable PCR-based detection and discrimination of BK and JC polyomaviruses. Souza LM, Savassi-Ribas F, Almeida SGS, Silva RNND, Baez CF, Zalis MG, Guimarães MAAM, Varella RB. Rev Inst Med Trop Sao Paulo. 2018;60:e47.
- 60 Simple technique for in field samples collection in the cases of skin rash illness and subsequent PCR detection of orthopox viruses and varicella zoster virus. Dumont C, Irenege LM, Magazani EK, Garin D, Muyembe JJ, Bentahir M, Gala JL. PLoS One. 2014 May 19;9(5):e96930

INVITEK DIAGNOSTICS GERMANY

Invitek Molecular GmbH
Robert-Roessle-Str. 10
13125 Berlin, Germany

Phone: +49 30 9489 2908

info@invitek.com

INVITEK DIAGNOSTICS PORTUGAL

ALS Life Sciences, SA
Zona Industrial de Tondela, ZIM II, Lote 6
3460-070 Tondela, Portugal

Phone: +351 232 817 817